

## NuclearIQ\_v1.3.0.3 Release Notes

### Issues Resolved within NuclearIQ

ID	Title
43761	Resolved the issue of the Method Definition not assigning a method to an analysis when the method has no Pre-QC requirement. These unassigned methods were not listed when the right click, Select Method was selected.
44070	Resolved issues when transferring gamma spectrum data to look only at the sample date/time, and allow the same system name and analysis date and time when from different unit numbers. The gamma spectrum data routine now looks at Site, Unit, and SystemName for a particular Sample date/time to ensure entries are unique.
44232	Resolved the issue of QC n-Number of points flag being applied to Non-Statistical Limits.
44285	Resolved the issue of the DBNull error received for each click on the Combination Plot Manager. Also, enabled the Method drop down list to be filtered by the selected site/unit/system/analysis.
44325	Resolved the issue on the Limit Evaluation Order form to correct the way the rows were highlighted when moved to new positions.
44327	Resolved the issue of the Method Definition screen showing a "cryptic" error message when attempting to delete a Test Stand, error message now states "Test Stand can not be deleted because it is associated to a Method on the Method Definition screen".
44328	Resolved the issue on the Input Test form to ensure comments associated with limit violations auto refresh on the form at save.
44330	Resolved the issue of the Rate of Change and Order of Magnitude inconsistently sending the appropriate warning message when exceeded.
44339	Resolved the issue of the Return to Spec email notifications not working when Notify on Warn is selected.
44347	Resolved the issue of the Corrective Action Guidelines (CAGs) not appearing on the Warning Limit notification screen.
44352	Resolved the issue of the Limits Effective Date not recognizing the time component.
44659	Resolved the issue of the error message received when more than one analysis is renamed on the Rename Analysis form.
44720	Resolved the issue of the error message received when a QC column is developed as a calculation column, and that column is the last column of the system.
44757	Resolved the issue of the application allowing QC data to be entered with future date/times.
44892	Resolved the issue of the error message received on the Method Definition form that was received after editing one method, to inform the user that another method is missing required information. This disables Save, but the user did not know what information was missing. This message was received even when no required information was missing.
44973	Resolved the issue of the Batch Scheduler having the ability to run twice within a single minute. Made the default timer 60 seconds to prevent this.
45284	Resolved the issue of the Evaluation Order form missing non-limit name entries.
45499	Resolved the issue of the incorrect position of the Input QC button on the Qc Input form. The button was renamed to Input New QC and positioned correctly.
45505	Resolved an issue with the Instrument Logbook truncating Calibration History entries when the Input Test form was open.

45506	Resolved an issue when using the right click, Enter/Review Analysis Comments. Comments were not viewable until manual reload of the input test form. Form now auto reloads to show the comment upon saving it.
46116	Resolved the issue on the Method Definition form to filter the Test Stand Name list by Site and Unit Number.
46146	Resolved an issue where an unnecessary error message is displayed when saving the Calibration History on the Instrument Logbook.
46715	Resolved an issue with the @Today variable when using the Batch Scheduler.
<b><i>New Features added to NuclearIQ</i></b>	
<b>ID</b>	<b>Title</b>
43870	Provided an enhancement to enable a user to filter and print the filtered grid from the Schedule Review form.
44315	Enhancements to the Instrument Logbook form. Test Stand Location was moved to the Test Stand Names form to allow the Location to be a drop down list on the Instrument Logbook form. Added filter boxes to the Log Book tab for Date/Author/Comment filters. Added filter boxes to the Calibration tab for Date/Login Name/Reason filtering.
<b><i>Updates made within NuclearIQ</i></b>	
<b>ID</b>	<b>Title</b>
44326	Resolved the issue of ordering Test Input List Items in a default ascending order.
44329	Resolved the issue of ordering Test Input List Item Library in a default ascending order for Group Names and then List Item Names.
44902	Resolved an issue with the Check Requirement Errors feature to check to see if the Requirement Mark date is valid.
44966	Updated the Batch Scheduler time adjustment spinners to limit months to 12, days to 31, hours to 23, and minutes to 59.
45496	Changed the default button focus on the schedule review form from the Close Schedule to the Close Form button.
45503	Enhanced the Plot Options to include Highlight Remarks (default True) and Tag Non Validated (default False).